

Installing Spectrum Viewer on Windows 7 & Windows 10

From the EDAX web site, **Support** tab and **Spectrum Viewer** link

Download the file

- 📁 If using Internet Explorer in Windows 7 or Edge in Windows 10, either right-click on the zip file link, or left click on it.
- 📁 If using a **Left-Click** on the file, a pop up with a prompt to Open or Save the file will be displayed. **Save** the file, first.
- 📁 Use your File Explorer to go to the Downloaded file directory.
- 📁 Then, **Extract all** files from the saved zip file on your local computer.

Left-click download

Service Support
Software Licensing
Spectrum Viewer
Technical Support
Training Schools

Contact EDAX

For Windows XP and older
[Spectrum Viewer Zip](#)
Modified: July 21, 2005
(481 KB)

For Windows 7 & 10
[Spectrum Viewer for Windows 7 & 10 Zip](#)
[Spectrum Viewer on Windows 7 Basic Instructions](#)
Modified: Feb. 2, 2017
(703 KB)

Spectrum Viewer 4.0 is a free utility that displays EDAX spectra (.spc files) on any PC, including multiple spectral overlays with X & Y offsets.

Follow EDAX on
LinkedIn Twitter WordPress YouTube

[Home](#) - [Careers](#) - [Contact Us](#)

© Copyright 2017 EDAX Inc. All Rights Reserved
[Privacy Policy](#) - [Trademarks](#) - [Terms & Conditions](#) - [Sitemap](#) - [www.ametek.com](#)

Spectrum_Viewer_Windows_7_10.zip finished downloading. Open Open folder View downloads X

Right-click download

If using a **Right-Click** on the file, choose **Save Target As....**

- 📁 Use your File Explorer to go to the Downloaded file directory.
- 📁 Then, **Extract all** files from the saved zip file on your local computer

Install the software

📁 Once the file is downloaded and files are extracted, you can run the install program from the Extracted files. Be sure to have **Administrator** rights.

📁 After extracting the files, right-click on **InstallToC.bat** and select to **Run as Administrator**.

📁 There should be three prompts showing a process succeeded:

```
C:\WINDOWS\System32\cmd.exe
QUIZ.SPC
ReadMe.txt
Sioxyn.spc
SpcViewer.exe
spectrum viewer installation on windows 7.pdf
Stnlss.spc
Views.dll
    9 file(s) copied.

C:\Users\khudson\Downloads\Spectrum_Viewer_Windows_7_10\Spectrum Viewer - Win7>c:\win
dows\system32\regsvr32.exe "c:\edax32\sys\Views.dll"

C:\Users\khudson\Downloads\Spectrum_Viewer_Windows_7_10\Spectrum Viewer - Win7>Rem Re
gister Display Component for Windows NT/

C:\Users\khudson\Downloads\Spectrum_Viewer_Windows_7_10\Spectrum Viewer - Win7>Rem Re
gister Display Component for Windows 95/98

C:\Users\khudson\Downloads\Spectrum_Viewer_Windows_7_10\Spectrum Viewer - Win7>C:\WIN
DOWS\System32\regsvr32 "c:\edax32\sys\Views.dll"
```

📁 The program will be installed in **c:\edax32\sys** folder.

📁 Run **SpcViewer** from the folder; **c:\edax32\sys**.

Do NOT run the SpcViewer from the Download or Extracted files folder.